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pH Calibration

For best results, periodic calibration with known, accurate standards is recommended. Calibrate with standards that bracket your intended measuring range while including a neutral standard (pH 7.00 or 6.86). For example, if you expect to measure samples from pH 6.2 to pH 9.5, calibration with 4.01, 7.00, and 10.01 standards will work well. Provide stirring for best results. After calibration with two or more points, the active slope segment of the measurement will be visible on the bottom display during measurement. 100 % slope will be shown if only one calibration point is performed and "- - -" if no calibration is performed. The meter will automatically return to measurement mode upon successful completion of the number of specified calibration points. To specify a different number of pH calibration points, See pH Buffer Options.

Using Automatic Buffer Recognition

1. While in pH measurement mode, dip the pH and ATC sensor(s) into your first standard, then press CAL. The primary display will search for the nearest standard value, while the secondary display will show the un-adjusted value.

2. When the "READY" indicator appears, press ENTER to accept. The primary reading will flash "DONE".

3. Rinse your electrode(s) then dip into the next pH standard. The primary display will search for the nearest standard value that has not yet been calibrated, while the secondary display will show the unadjusted value. When the "READY" indicator appears, press ENTER to accept.

4. To calibrate another pH standard repeat Step 3 or press MEAS to return to pH measurement mode.