**Bedienungsanleitung**  
**Operating Instructions**  
**Mode d'emploi**

- » Star-Thermo Exsikkator
- » Star-Thermo Desiccators
- » Dessiccateur Thermo-Star

### SICCO Star-Thermo-Exsikkatoren, Artikelnummer: V 1830-07

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### SICCO Star-Thermo Desiccators, Cat.-No.: V 1830-07

#### OPERATING INSTRUCTIONS

1. Supplied with	8
2. Security Advice	8
3. Taking into operation	8
4. Mode of function	11
5. Service / Maintenance	11
6. Disturbances / Repairs	11
7. Waste Disposal	12
8. Specifications	12
9. Accessories	12

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Star-Thermo Desiccators are made for storing your samples and products at a constant temperature. The special 2-panes insulation (inside glass, outside PMMA) reduces heat emission and the solid aluminium frame provides maximum security.

The Star-Thermo Desiccators are ideal for warming your samples and substances in order to reduce the reduction times or for temporary storage at a constant ambient temperature.

# Sicco Star-Thermo Desiccators

## 1. Supplied with

- 1 SICCO Star-Thermo Desiccator
- 4 Shelves
- 8 Rails for Shelves
- 1 Operating Instructions

## 2. Security Advice

- » Do not take a damaged desiccator into operation.
- » Do not put any explosive or inflammable products into the desiccator.
- » Do not store any solvents or acids in the desiccator.
- » Do not store any explosive or inflammable products near the desiccator. Keep a distance of at least one meter.
- » During the operation the desiccator will get hot. ATTENTION: Risk of burns.
- » SICCO Star-Thermo Desiccators are not suitable for overpressure or vacuum.
- » Provide a safe stand for the desiccator.
- » SICCO Star-Thermo Desiccators are to be used in closed and dry rooms only.
- » Protect the Star-Thermo Desiccator from splash water.
- » Place the desiccator in a reachable working height.
- » The ambient temperature range of  $-20\text{ }^{\circ}\text{C}$  to  $+70\text{ }^{\circ}\text{C}$  shall not be exceeded.
- » Do not expose the desiccator to direct sunlight.
- » Remove only one shelf at once.
- » The temperature inside the desiccator shall not exceed  $+60\text{ }^{\circ}\text{C}$ .
- » Do not put any products into the desiccator which will fly about due to the air circulation, e.g. powdery substances or light membranes.
- » Replace the microfuse only with an equal 4 ampère fuse.

## 3. Taking into operation

- » Remove the protection cover.
- 
- » Remove the protection packing of the power plug. Do not use a knife. Plug the power plug in a 230 V outlet.


## Sicco Star-Thermo Desiccators

» Remove the protection film from the shelves.


» Opening the door: Hold the desiccator housing with one hand while you open the door by pulling the door grip with the other hand.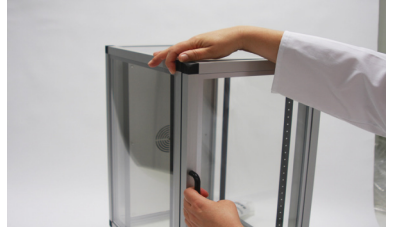


» **Please note:** Due to the strong magnetic lock, it is harder to open the door the longer it was closed before. This is normal and can be compensated by pulling stronger than in the everyday use.

» The temperature sensor (1) can be positioned freely by attaching it to the suspension arrangement (2). That way you can monitor the temperature at the required point.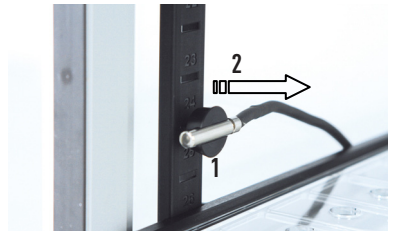


» To change the position, remove the retainer with the temperature sensor (1) from the suspension arrangement (2).


» Pull the cable (1) of the temperature sensor carefully upwards in a vertical direction.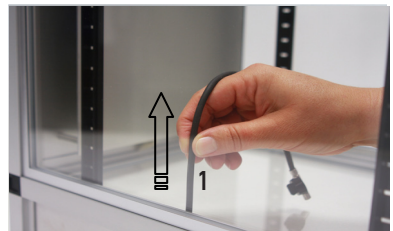

» Once the cable is long enough to reach the required position you can attach the temperature sensor (1) by clicking it into the suspension arrangement (2).

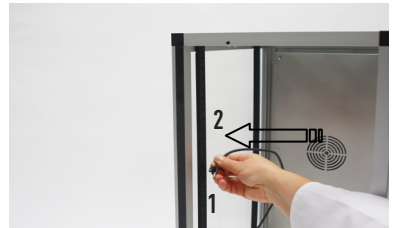
» Mount the rails (3) horizontally in pairs on both sides of the suspension arrangement (2). The engagement of the rail (3) is clearly audible. For a better orientation, the levels are marked by numbers on the suspension arrangement (2).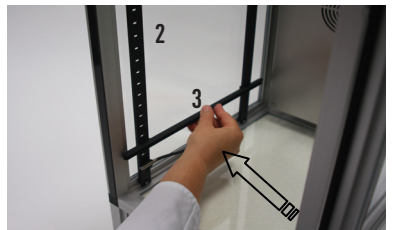


## Sicco Star-Thermo Desiccators


- » The shelves (4) can then be slid in on the rails (3).
- » Mount another pair of rails (3) and slide in another shelf (4).
- » **Please note:** Mount the rails in pairs and slide in immediately a shelf to avoid that other rails can fall down accidentally.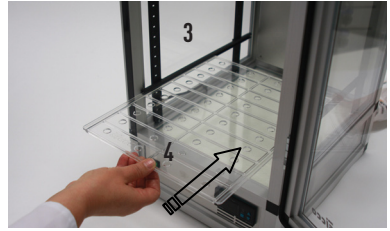

- » Close the door.
- » Turn on the power switch (5) of the desiccator. The display shows the indoor temperature.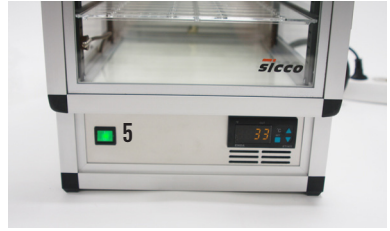

- » **Attention:** To reduce a drop of temperature, a contact switch (6) inside the door frame turns off automatically the heating and the fan when the door is open. If you close the door the desiccator will heat again with the set value.

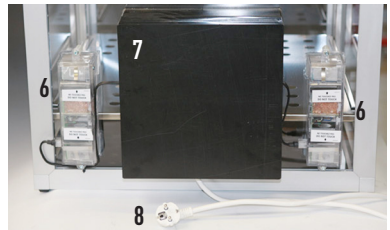- » For adjusting the temperature press the "SET" key permanently. The display shows the set temperature. By pressing simultaneously the "SET" key and one of the "arrow" keys you can change the set temperature.
- » For exact regulation the set temperature has to be 5°C above room temperature.
- » During operation the set temperature can be displayed by pressing the "SET" key. The temperature display flashes.


## 4. Mode of function

The fan on the back side sucks in the air inside the desiccator and blows it over the heating spiral. At the same time, the warmed air flows back into the desiccator. This air circulation provides an even heating of the interior space.

The temperature sensor inside monitors continuously the actual temperature. Once the temperature has reached the set value, the heating is switched off. The fan keeps on working for providing best heat distribution.

As soon as the internal temperature drops 2°C below the set value, the heating is switched on again automatically.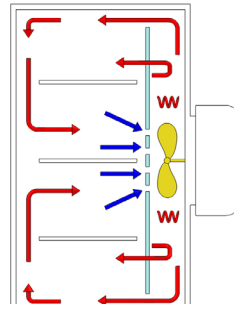

## 5. Service / Maintenance

- » Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.
- » Keep the magnetic door strip clean.

## 6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then, a solution to your problem cannot be found, we will ask you to return the cleaned desiccator in the original packaging to us. Then we will check whether a cost-effective repair is possible.

Objection	Reason and help
Die eingestellte Soll-Temperatur wird nicht erreicht.	<ul style="list-style-type: none"> <li>» The door is not closed tightly. The heating and the fan are turned off automatically by the contact switch(6) inside the door frame. Close the door.</li> <li>» The set temperature is lower than the current room temperature. Make sure that the set temperature is at least 5 °C above the current room temperature.</li> </ul>
No reaction when the power switch is turned on.	<ul style="list-style-type: none"> <li>» The power plug is not plugged in. Plug it in a 230 V outlet.</li> <li>» The microfuse is broken. Replace the microfuse as Follows.</li> </ul>


### How to replace the microfuse:

- » Unplug the power plug from the outlet.
- » Open the fuse housing (7). Use a screwdriver with a width of 5,5 mm and screw off the housing counterclockwise.

## Sicco Star-Thermo Desiccators

- » Replace the microfuse (8) with an equal 4 ampère fuse.  
Screw in the fuse housing (7) with the new microfuse clockwise.
- » **Attention:** In case that the new microfuse is broken after turning on the desiccator again, please contact an electrician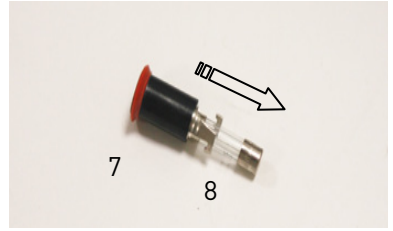

### 7. Waste Disposal

The desiccators as well as their packing are made of recyclable materials. Please take further information about the materials used from the following specifications.

### 8. Specifications

#### SICCO Star-Thermo Desiccator

Cat.-No.:	V 1830-07
Material panels:	Inside glass, outside PMMA (Polymethylmethacrylat)
Temperature resistance:	- 20° bis + 70°C
Material frame:	Aluminium with connectors made of polyamide
Material of rails:	Polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D)	310 x 630 x 420 mm, additional place requirement for heating: 80 mm back side
Usable interior space (W x H x D):	260 x 480 x 330 mm
Weight:	approx. 16 kg
Maximum number of shelves:	26
Maximum all-over load per shelf	2,5 kg
Maximum all-over load:	10 kg
Temperature setting:	Room temperature plus 5 °C up to max. + 60 °C
Setting accuracy:	1 °C
Deviation of temperature:	approx. +/- 1,75 °C at + 40°C
Heating system:	air circulation
Power consumption:	850 W
Power connection:	230 V / 50 Hz
Fuse:	4 A microfuse

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BOHLENDER GmbH